

MST CUTICLE BLUE

USAGE DIRECTIONS:

(Wear rubber gloves to avoid staining hands. Use clean suitable container to test eggs. Return fluid to original container / seal for next use)

ACCIDENTAL SPILLAGE: absorb with sponge etc, finish rinse with clean water.

1. Submerge egg(s) for 1 minute
2. Remove eggs and rinse under running water (2 or 3 seconds) to remove surplus
3. Allow eggs to dry / monitor results

Results:

Full colour	= Good cuticle
Random colouring	= Irregular cuticle
Damage lines	= Handling problems
No colour	= No cuticle

Additional data information:

Internal membrane stain= Poor shell and or cuticle structure
(Use recommended chemicals when egg sanitising)

INGREDIANTS: NON HAZARDOUS

Blended mixture of cuticle sensitive dyes suspended in aqueous solution (or shipped in POWDER form) for non destructive egg testing.

Expiry Date: END FEB 2010

Batch No : MS 060208

MS TECHNOLOGIES LTD

Telford Way, Kettering, Northants, NN16 8UN, England

Tel: +44 (0) 1536 516778 Fax: +44 (0) 1536 310027

MS TECHNOLOGIES (USA) PIKEVILLE, TN 37367 Ph: (423) 881 3880

Shipping Method: FLUID FORM : READY TO USE
POWDER FORM : (ADD POTABLE WATER)
* ADD POTABLE WATER TO VOLUME
** TRANSFER POWDER CONTENT TO SUITABLE SIZE CONTAINER AND ADD POTABLE WATER TO VOLUME

CONTENTS:

1 LTR * 5 LTR * 25 LTR **